



Owner: [REDACTED]	Number: Bio-BPC-Form-11
Author: [REDACTED]	Version: 1
Active: 15/06/2021	Review: <QPulse_DocReviewDate>
Title: Appendix 1 - Fragment Analyzer Worksheet - Pfizer COMIRNATY	

Worksheet for Fragment Analyzer - RNA Integrity

Test Details			
SOP QPulse #	Bio-BPC-Method-26	Analysist	[REDACTED]
TRIM link to data files	D21-2862801, D21-2862802	Test Date	20/07/2021

Pipettes & Equipment	
Name	LIMS#
30-300 µL 12 channel pipette	-
p10 pipette	-
p50 pipette	-
p100 pipette	32792
p200 pipette	32892
Thermomixer	32660
Thermocycler	32865
P20	32891

Reagents & Consumables			
Details	Catalog #	Lot/Batch Number	Expiry date
48-Capillary Array, short 33 cm	A2300-4850-3355	022621-27sfs	-
Standard Sensitivity (SS) RNA kit Part 1 stored at 2-8°C	DNF-471-0500	6591350	8/12/2021
<i>Extra Blank solution</i>	<i>DNF-300-0008</i>	-	-
Standard Sensitivity (SS) RNA kit Part 2 stored at -20°C (Diluent Marker & Intercalating dye)	-	6595832	8/03/2022
<i>Extra Diluent marker</i>	<i>DNF-369-0004</i>	-	-
Standard Sensitivity (SS) RNA kit Part 3 stored at -70°C (RNA Ladder)	DNF-382-U020	6602189	6/04/2022
Capillary conditioning solution	DNF-475-0100	6588180	29/01/2022
DEPC water	AM9961	2004017	-
20% Triton-X100 / 30% Ethanol solution	In house	MC1JUN21-01	1/09/2021
RNA Separating Gel	DNF-265-0500	6597220	15/04/2022
Enter text.	Enter text.	Enter text.	Enter a date.
Enter text.	Enter text.	Enter text.	Enter a date.

Reagent Preparation			
REAGENT	STORAGE	Date Prepared	Expiry
Inlet Buffer - 10 mL 5X inlet buffer (fridge) - 40 mL MilliQ Water	Deep well plate, RT (Drawer B)	20/07/2021	21/07/2021 Prepare fresh
Rinse Buffer 200 µL per well of 0.25X TE buffer (stored in fridge, allow to come to RT) Can also use DEPC water	96 well plate, RT (Drawer M)	20/07/2021	21/07/2021 Prepare fresh
Capillary Storage Buffer 100 µL per well capillary storage solution (RT)	96 well plate, RT (Drawer 3)	16/07/2021	30/07/2021 2 weeks
Capillary Conditioning solution - 6 mL 5X capillary conditioning Soln (RT) - 24 mL Milli-Q Water	250 mL tube, RT (Capillary conditioning line)	20/07/2021	3/08/2021 2 weeks
RNA Separation gel - 23 mL RNA separating gel (fridge) - 2.3 µL Intercalating dye (-20°C)	50 mL tube, RT (Gel line 2)	20/07/2021	22/07/2021 48 hours
Empty waste tray and waste bottle Reagents can be scaled up if required – this table provides the minimum for a single run. Keep all samples, RNA ladder (-70°C), and RNA diluent marker (-20°C) on ice until ready for use Allow separating gel, blank solution, rinse buffer, inlet buffer (all stored in fridge), and the intercalating dye (-20°C) to come to RT before use Blank solution can be replaced with diluent marker. All wells need to contain a peak for capillary alignment.			

96 well Plate Layout

	1	2	3	4	5	6	7	8	9	10	11	12
A	BF-25	BF-25	BF-25	BF-25	BF-25	S6a	S5a	S4a	S3a	S2a	S1a	RM
B	BF-25	BF-25	BF-25	BF-25	BF-25	S6b	S5b	S4b	S3b	S2b	S1b	RM
C	BF-25	BF-25	BF-25	BF-25	BF-25	S6c	S5c	S4c	S3c	S2c	S1c	RM
D	BF-25	BF-25	BF-25	BF-25	BF-25	BF-25	BF-25	BF-25	BF-25	BF-25	BF-25	L

S1-6 = Samples in triplicate (a, b or c),
RM = Reference material
BF-25 = Blank solution provided in kit,
L = RNA ladder

This worksheet assumes the maximum of 6 samples per test. Any samples not included in the test must be crossed off the plate layout, and results table below

System Suitability Criteria – RNA Ladder			
Plate location (wells)	Choose an item.		
Parameter	Limits	Results	Comments
RNA ladder profile	Visually comparable to figure 4 of SOP	ok	PASS
All peaks present	15 200 500 1000 1500 2000 3000 4000 6000 nt	ok	PASS
Peak heights	<60000 RFU	ok	PASS
Assay Acceptance Criteria – Reference Material			
Plate location (wells)	A12 B12 C12		
LIMS #	2103001141		
BATCH #	EE8493		
EXPIRY	5/08/2021		
Parameter	Limits	Results	Comments
Profile	Visually comparable to DP electropherogram in SOP	Ok/ok/ok	PASS
Migration time	Approximately comparable to profile in SOP	4097/4097/4097	PASS
Lower marker present	LM peak	Ok/ok/ok	PASS
Peak heights	5000-600000 for 2/3 replicates	13860/12940/12365	PASS
No negative peaks or baseline shifts	No significant peaks/shifts	Ok/ok/ok	PASS
Reference Material Dilutions / Calculation / Notes			
thaw date: 20/07/21 270ng/uL = 20 uL of 530 ng/uL master stock + 19 uL DEPC water 90 ng/uL = 20 uL of 270 ng/uL + 40 uL 20%Tx100 solution			

Sample 1 Details	
Plate location (wells)	A11 B11 C11
LIMS #	2107002694
BATCH #	FE3430
EXPIRY	31/10/2021

Sample Acceptance Criteria			
Parameter	Limits	Results	Comments
Migration time	Comparable to RM	4097/4097/4097	PASS
Lower marker	LM must be present	Ok/ok/ok	PASS
Peak heights	5000-60000 RFU for 2/3 replicates	15646/17433/17849	PASS

Test Results					
Parameters	Limits	Results			Comments
		Average	SD	%RSD	
% RNA Integrity		67.7	1.66	2.46	PASS
% Late Migrating Species		2.97	0.15	5.15	

Sample Dilutions / Calculation / Notes
Thaw date – 19/07/2021 – stored in cold room 270ng/uL = 20 uL of 500 ng/uL master stock + 17uL DEPC water 90 ng/uL = 20 uL of 270 ng/uL + 40 uL 20%Tx100 solution

Sample Results	
PASS	
Analysist	
Checked by	Enter text.

Sample 2 Details	
Plate location (wells)	Choose an item.
LIMS #	Click or tap here to enter text.
BATCH #	Click or tap here to enter text.
EXPIRY	Enter date.

Sample Acceptance Criteria			
Parameter	Limits	Results	Comments
Migration time	Enter text.	Enter text.	Choose an item.
Lower marker	Enter text.	Enter text.	Choose an item.
Peak heights	Enter text.	Enter text.	Choose an item.

Test Results					
Parameters	Limits	Results			Comments
		Average	SD	%RSD	
% RNA Integrity	Enter text.	Enter text.	Enter text.	Enter text.	Choose an item.
% Late Migrating Species		Enter text.	Enter text.	Enter text.	

Sample Dilutions / Calculation / Notes
Enter text.

Sample Results	
Choose an item.	
Analysist	Enter text.
Checked by	Enter text.

Sample 3 Details	
Plate location (wells)	Choose an item.
LIMS #	Click or tap here to enter text.
BATCH #	Click or tap here to enter text.
EXPIRY	Enter date.

Sample Acceptance Criteria			
Parameter	Limits	Results	Comments
Migration time	Enter text.	Enter text.	Choose an item.
Lower marker	Enter text.	Enter text.	Choose an item.
Peak heights	Enter text.	Enter text.	Choose an item.

Test Results					
Parameters	Limits	Results			Comments
		Average	SD	%RSD	
% RNA Integrity	Enter text.	Enter text.	Enter text.	Enter text.	Choose an item.
% Late Migrating Species		Enter text.	Enter text.	Enter text.	

Sample Dilutions / Calculation / Notes
Enter text.

Sample Results	
Choose an item.	
Analysist	Enter text.
Checked by	Enter text.

Sample 4 Details	
Plate location (wells)	Choose an item.
LIMS #	Click or tap here to enter text.
BATCH #	Click or tap here to enter text.
EXPIRY	Enter date.

Sample Acceptance Criteria			
Parameter	Limits	Results	Comments
Migration time	Enter text.	Enter text.	Choose an item.
Lower marker	Enter text.	Enter text.	Choose an item.
Peak heights	Enter text.	Enter text.	Choose an item.

Test Results					
Parameters	Limits	Results			Comments
		Average	SD	%RSD	
% RNA Integrity	Enter text.	Enter text.	Enter text.	Enter text.	Choose an item.
% Late Migrating Species		Enter text.	Enter text.	Enter text.	

Sample Dilutions / Calculation / Notes
Enter text.

Sample Results	
Choose an item.	
Analysist	Enter text.
Checked by	Enter text.

Sample 5 Details	
Plate location (wells)	Choose an item.
LIMS #	Click or tap here to enter text.
BATCH #	Click or tap here to enter text.
EXPIRY	Enter date.

Acceptance Criteria			
Parameter	Limits	Results	Comments
Migration time	Enter text.	Enter text.	Choose an item.
Lower marker	Enter text.	Enter text.	Choose an item.
Peak heights	Enter text.	Enter text.	Choose an item.

Test Results					
Parameters	Limits	Results			Comments
		Average	SD	%RSD	
% RNA Integrity	Enter text.	Enter text.	Enter text.	Enter text.	Choose an item.
% Late Migrating Species		Enter text.	Enter text.	Enter text.	

Sample Dilutions / Calculation / Notes
Enter text.

Sample Results	
Choose an item.	
Analysist	Enter text.
Checked by	Enter text.

Sample 6 Details	
Plate location (wells)	Choose an item.
LIMS #	Click or tap here to enter text.
BATCH #	Click or tap here to enter text.
EXPIRY	Enter date.

Acceptance Criteria			
Parameter	Limits	Results	Comments
Migration time	Enter text.	Enter text.	Choose an item.
Lower marker	Enter text.	Enter text.	Choose an item.
Peak heights	Enter text.	Enter text.	Choose an item.

Test Results					
Parameters	Limits	Results			Comments
		Average	SD	%RSD	
% RNA Integrity	Enter text.	Enter text.	Enter text.	Enter text.	Choose an item.
% Late Migrating Species		Enter text.	Enter text.	Enter text.	

Sample Dilutions / Calculation / Notes
Enter text.

Sample Results	
Choose an item.	
Analysist	Enter text.
Checked by	Enter text.

Notes

Enter text.